

## **Activate or Register QCapture Pro 7™ on a Computer without Internet Access**

Verify the computer is running one of the operating systems recommended for QCapture Pro 7 Software:

- Windows® XP SP 3 and later (32-bit)
- Windows® Vista Business and Ultimate SP 2 and later (32 and 64 bit)
- Windows® 7 Professional and Ultimate (32 and 64 bit)

You will need the following:

- QCapture Pro 7 already installed and running the 30 day trial
- QCapture Pro 7 USB key
- A USB flash drive

### **Instructions**

1. Insert USB flash drive
2. Insert USB key for QCapture Pro 7
3. Start QCapture Pro 7
4. You will see the following screen (Figure 1), click *OK*

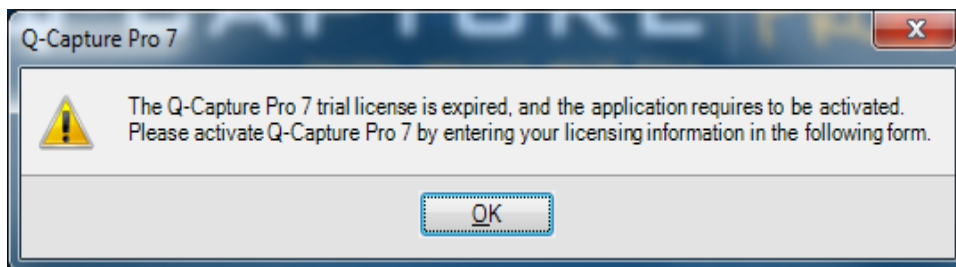


Figure 1

19535 56<sup>th</sup> Avenue, Suite 101

Surrey, BC

Canada V3S 6K3

tel 604.530.5800

fax 604.648.8277

[www.qimaging.com](http://www.qimaging.com)

5. Click *Cancel* on the next window (Figure 2).  
*Note the Error as the software is trying to access the internet to check for updates on your PC that is currently not network connected.*

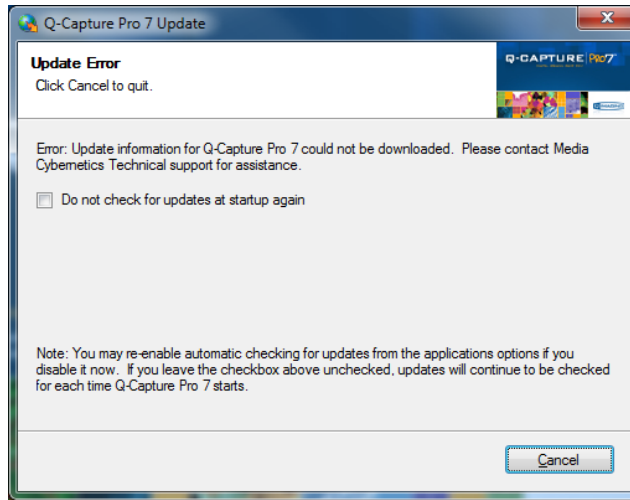


Figure 2

6. Choose a USB drive and click *OK* (Figure 3) and proceed to click *OK* (Figure 4).

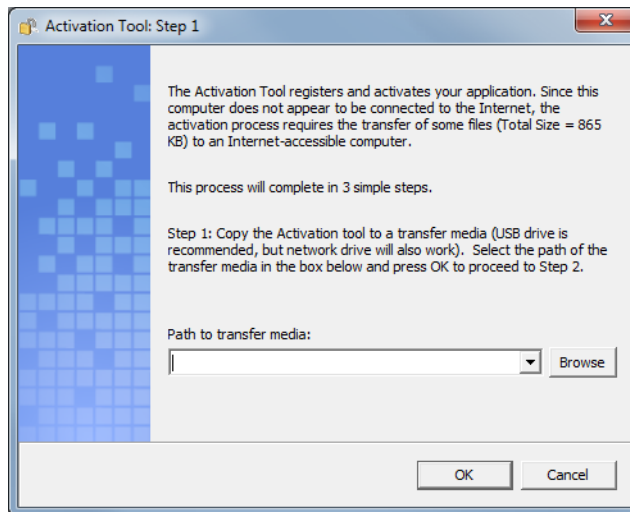


Figure 3

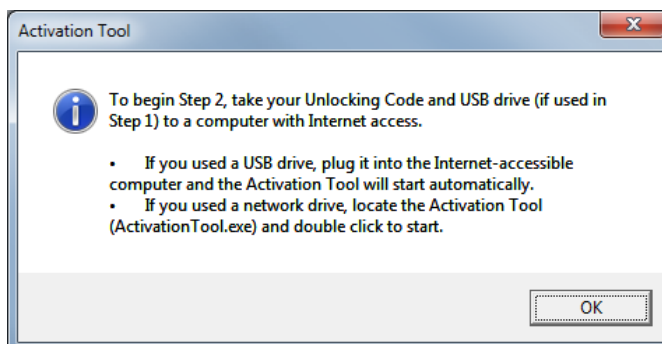


Figure 4

7. Install the USB flash drive on a computer with internet access.
8. Browse the USB flash drive on the computer with internet access and double click on *ActivationTool* (Figure 5) unless it already started by itself.

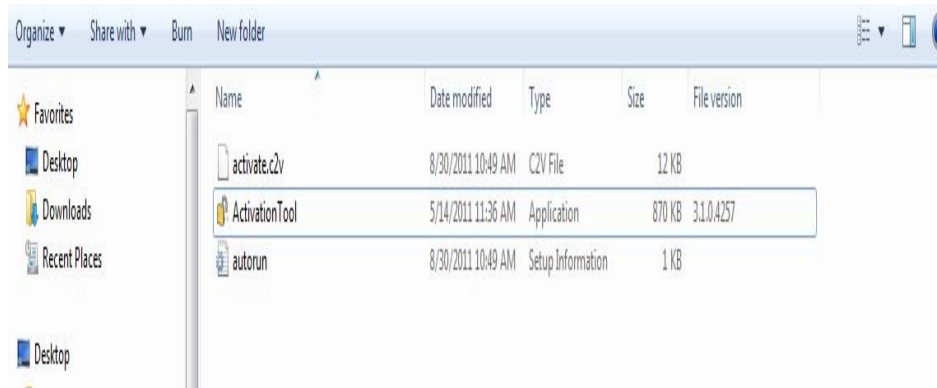


Figure 5

9. Complete the form and submit (Figure 6), then click *OK* (Figure 7).

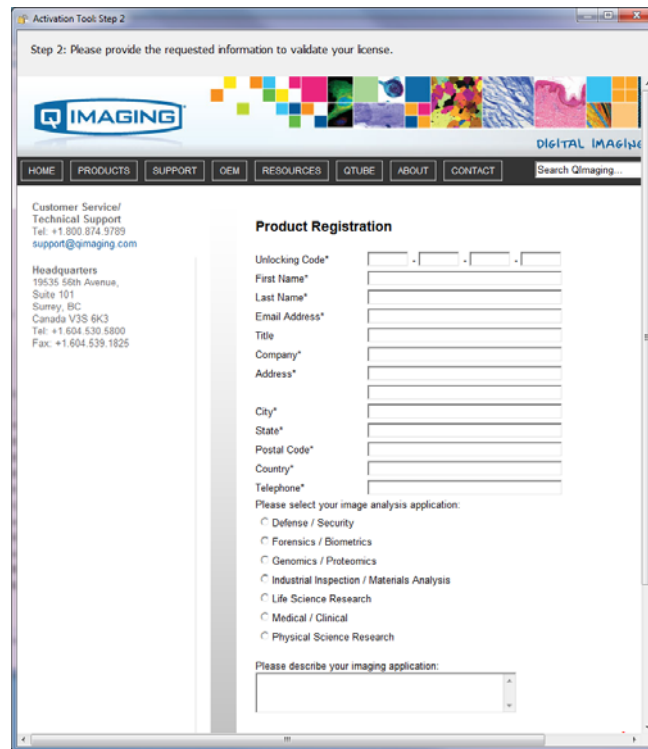


Figure 6

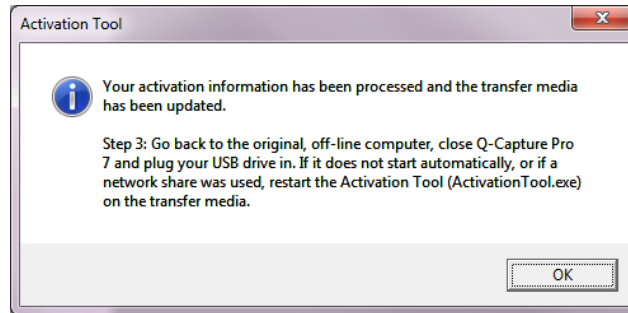


Figure 7

10. Click *OK*. Now the USB flash drive will have one extra file: *Q-Capture Pro 7.isn* (Figure 8)

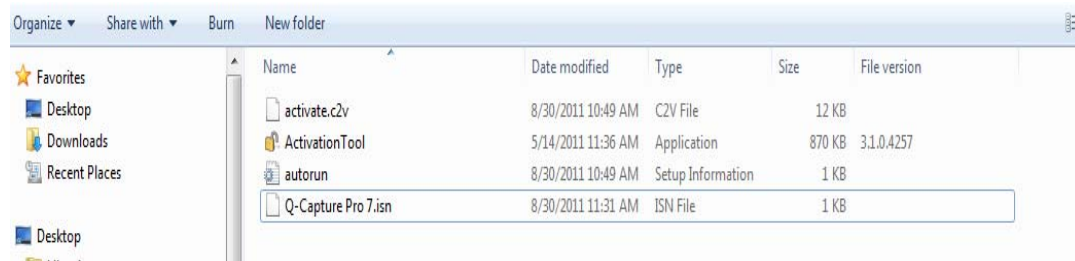


Figure 8

11. Remove the USB flash drive and install it on the computer without internet access.
12. If the activation does not start, run the *ActivationTool* on the USB key. You will see the following message (Figure 9).

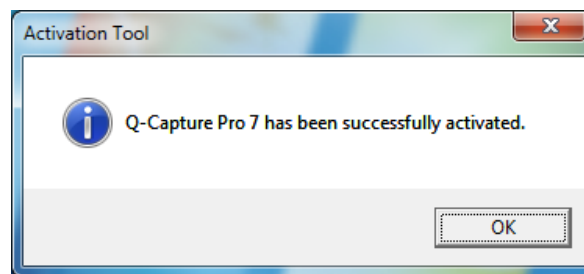


Figure 9

13. Click *OK*
14. QCapture Pro 7 is now activated.